

ACR17 Drawer Instructions

This pharmacy refrigerator can hold up to nine drawers, but it is best to use no more than seven to make sure air is evenly distributed throughout the interior. You can mix shelves and drawers as needed.

How to Install the ACR17 Drawer

1. Each drawer ships with two **interior dividers** for optional storage organization. If using these dividers, install them inside the drawer **BEFORE** inserting the drawer into the refrigerator. To install, place the divider diagonally inside the drawer, hooking one side in the desired position. Slide the other side of the divider in until it snaps into place.



Note: If drawers are being installed towards the bottom of the refrigerator, start at the fourth notch up.

2. **Align** the hooks of the bracket at a 45° angle upwards and push down to secure in place. You can use the number or letter code printed above the holes to make sure the brackets are lined up evenly.



3. **Slide** the drawer into the tracks one side at a time by guiding the drawer into the track until you hear a click. Repeat this same step for the other side. Slide the drawer in and out to make sure it is correctly aligned and leveled.

NOTE: If the drawer gets stuck or does not feel level, **DO NOT** try to push it in, as you may damage the glides. Instead, remove the drawer (see next page for instructions) and try again, slowly aligning each side and gently pushing the drawer into place.



Removing the ACR17 Drawer

1. Empty all stored contents from the drawer before removing it from the unit.
2. Pull the drawer all the way out.
3. You will see a small black latch on each side of the drawer. Push the latch on the left side up and pull the latch on the right side down at the same time. Proceed to slide the drawer out.
4. To remove the tracks, simply tilt each bracket up and pull the track out.



LEFT SIDE



RIGHT SIDE