3-DRAWER ALL-FREEZER
for Medical and General-Purpose Storage

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Use & Care Guide

Mode d'emploi

BEFORE USE, PLEASE READ AND FOLLOW
ALL SAFETY RULES AND OPERATING INSTRUCTIONS

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FELIX STORCH, INC.
An ISO 9001:2015 registered company
770 Garrison Avenue
Bronx, New York 10474
www.summitappliance.com

Table of Contents

Important Safeguards	2 - 3
Installation	4 - 5
<i>General Instructions</i>	4
<i>Location and Clearances</i>	4 - 5
<i>Electrical Connection</i>	5
Operation	6 - 7
<i>When Using the Appliance</i>	6
<i>Removal/Replacement of the Drawers</i>	6
<i>Temperature Selection and Monitoring</i>	7
<i>Defrosting</i>	7
Appliance Care	8 - 9
<i>Cleaning the Appliance</i>	8
<i>Loss of Power</i>	8
<i>Moving the Appliance</i>	9
<i>Disposing of a Worn-Out Appliance</i>	9
Troubleshooting	10
Limited Warranty	11
French version of this manual	13 - 23

IMPORTANT SAFEGUARDS

Your safety and the safety of others are very important to us.

Before the appliance is used, it must be properly positioned and installed as described in this manual, so read the manual carefully. To reduce the risk of fire, electrical shock or injury when using the appliance, follow basic precautions, including the following:

DANGER

- Plug into a grounded 3-prong outlet, do not remove grounding prong, do not use an adapter, and do not use an extension cord.
- Replace all panels before operating.
- It is recommended that a separate circuit serving only your appliance be provided. Use receptacles that cannot be turned off by a switch or pull chain.
- Never clean appliance parts with flammable fluids. These fumes can create a fire hazard or explosion. And do not store or use gasoline or other flammable vapors and liquids in the vicinity of this or any other appliance. The fumes can create a fire hazard or explosion.
- Do not connect or disconnect the electric plug with wet hands.
- Unplug the appliance or disconnect power before cleaning or servicing. Failure to do so can result in electrical shock or death.
- Do not attempt to repair or replace any part of your appliance unless it is specifically recommended in this manual. All other servicing should be referred to a qualified technician.

WARNING

FOLLOW WARNING CALLOUTS BELOW ONLY WHEN APPLICABLE TO YOUR MODEL

- Use two or more people to move and install appliance. Failure to do so can result in back or other injury.
- To ensure proper ventilation for your appliance, the front of the unit must be completely unobstructed. Choose a well-ventilated area with temperatures above 44°F (7°C) and below 95°F (35°C).
- The appliance should not be located next to ovens, grills or other sources of high heat.
- The appliance must be installed with all electrical, water and drain connections in accordance with state and local codes. A standard electrical supply (115 V AC only, 60 Hz), properly grounded in accordance with the National Electrical Code and local codes and ordinances, is required.
- Do not kink or pinch the power supply cord of the appliance.

- The size of the fuse (or circuit breaker) should be 15 amperes.
- It is important that the appliance be leveled in order to work properly. You may need to make several adjustments to level it.
- All installations must be in accordance with local plumbing code requirements.
- Make certain that the pipes are not pinched, kinked or damaged during installation.
- Check for leaks after connection.
- Never allow children to operate, play with or crawl inside the appliance.
- Do not use solvent-based cleaning agents or abrasives on the interior. These cleaners may damage or discolor the interior.
- Use this appliance only for its intended purpose as described in this *Use & Care Guide*.
- Keep fingers out of the “pinch point” areas. Clearances between the drawers and cabinet are necessarily small. Be careful closing drawers when children are in the area.

Risk of child entrapment!

Child entrapment and suffocation are not problems of the past. Junked or abandoned appliances are still dangerous, even if they will “just sit in the garage for a few days”.

Before discarding your old appliance:

- *Take off the doors*
- *Leave the shelves in place so that children may not easily climb inside.*

- SAVE THESE INSTRUCTIONS -

Installation

Excessive Weight Hazard !

***Use two or more people to move and install your appliance.
Failure to do so can result in back or other injury.***

General Instructions:

1. Remove packaging materials.

2. **CAUTION:**

*After unpacking you **MUST** allow this freezer to stand upright for at least 2 hours to allow the lubricant and refrigerant to drain back into the compressor and stabilize. Failure to do so may adversely affect performance and the lifetime of this unit.*

3. Remove tape and glue residue from surfaces before turning on the appliance. Rub a small amount of liquid dish soap over the adhesive with your fingers. Rinse with warm water and dry.

4. Do not use sharp instruments, rubbing alcohol, flammable fluids, or abrasive cleaners to remove tape or glue. These products can damage the surface of your appliance.

5. **When Moving Your Appliance:**

Your appliance is heavy. When moving the appliance for cleaning or service, be sure to protect the floor. Always pull the appliance straight out when moving it. Do not wiggle or “walk” the appliance when trying to move it, because floor damage can occur.

6. **Clean Before Using:**

After you remove all of the packaging materials, clean the inside of your appliance before using it. See cleaning instructions in *Appliance Care*.

7. Some models have leveling legs. Level as necessary for best operation.

8. If your appliance has casters, after moving into place, lock the brakes on those casters supplied with brakes. Always unlock before moving the appliance for cleaning.

Location and Clearances:

IMPORTANT: Do not install the appliance near an oven, radiator, or other heat source, nor in a location where the temperature will fall below 41°F (5°C). Do not place in direct sunlight.

Check the strength of the floor before installing your appliance. It should hold

the weight of a fully loaded appliance.

To ensure optimum operation and proper ventilation, for free-standing installation, allow at least 4" (10 cm) of airspace at the top and on both sides of the appliance and also 4" between the back of the appliance and the wall. If the freezer is to be built-in, allow 3/16" clearance on both sides, 2" at the rear and 1/16" at the top. Be sure that airflow under the appliance is not blocked.

If your appliance was supplied with anti-tip brackets, be sure to install them following the directions included with the brackets.

Electrical Connection:

A 115 Volt (AC only) 60 Hz grounded electrical supply with a 15- or 20- amp fuse or circuit breaker is required. Have the wall outlet and circuit checked by a qualified electrician to make sure the outlet is properly grounded. When a standard 2-prong wall outlet is encountered, it is your responsibility and obligation to have it replaced with a properly grounded 3-prong wall outlet.

It is recommended that a separate circuit serving only your appliance be provided.

Use outlets that cannot be turned off by a switch or pull chain.

DO NOT USE AN EXTENSION CORD.

Electrical Shock Hazard

- ***Plug into a grounded 3-prong outlet.***
- ***Do not remove ground prong.***
- ***Do not use an adapter.***
- ***Do not use an extension cord.***

Failure to follow these instructions can result in fire, electrical shock or death.

Operation

When Using the Appliance:

In order to ensure proper temperature, allow proper airflow in and around the appliance. When stocking the freezer, do not block off any air vents. If the vents are blocked, airflow will be hindered, and temperature and moisture problems may occur. See page 5 for proper clearances around the unit.

Let the freezer cool for at least 6 to 8 hours before adding contents. Allow 24 hours for the temperature to stabilize.

NOTE: When the appliance is first started, the compressor will run constantly until the appliance is cooled. It may take 6 to 8 hours (or longer), depending on room temperature and the number of times the appliance drawers are opened. After the freezer has cooled, the compressor will turn on and off as needed to keep the appliance at the proper temperature.

For optimum temperature performance, it is recommended that this unit be at least 30% full during use. If you are only using this unit to store small loads, we recommend adding some freezer packs to ensure temperature stability.

The outside of the freezer may feel warm. This is normal. The appliance's design and main function is to remove heat from packages and air space inside the appliance. The heat is transferred to the room air, making the outside of the appliance feel warm. *(In some models the condenser that gets warm is just inside the side walls, making these walls warm to the touch.)*

ENERGY-SAVING TIPS

- The more airflow provided around an appliance, the less energy it will use.
- Keep ice build-up inside the freezer to less than 1/2". At this point you should manually defrost the unit. See page 7.
- Keep drawers open a minimum of time, and shut securely.

Removal/Replacement of the Drawers:

To remove a drawer from the freezer:

- Pull the drawer out towards you completely.
- On both sides of the drawer there are hinges.
- On each hinge there is a black lever.
- The left lever should be pulled up.
- The right lever should be pulled down.
- Both levers should be pulled at the same time while continuing to pull out the drawer.

To reinstall the drawers:

- Line up the hinges on the drawer with the hinges on the appliance and gently push the drawer in.
- Once it is lined up, close the drawer completely then open again to make sure the drawer is on track.

Temperature Selection and Monitoring:

The temperature inside the freezer is controlled by a digital thermostat located at the bottom of the unit at the right end of the front grille. The LED display normally shows the temperature inside the freezer. If you find it necessary to change the temperature setting (*Set Point*) of your appliance, refer to the following instructions. The front panel of the thermostat is illustrated here:

To see the Set Point:

Push and immediately release the *SET* key. The display will show the *Set Point* value. Wait 5 seconds and the interior temperature value will display again.

To change the Set Point:

Push and hold the *SET* key for more than 3 seconds to change the *Set Point* value. The value of the *set point* will be displayed, and the “snowflake” LED will start blinking.

To change the SET value:

Push the UP or DOWN arrow to the desired setting. The new setting will be memorized after 15 seconds, provided no other keys are pushed.

NOTE: The minimum and maximum range of your thermostat is preset and locked by us so it cannot easily be accidentally changed. To reset this range lock, either call us or find information on our website under *Troubleshooting*.

Defrosting:

The EQFM3D series are manual defrost all-freezers. No internal fans are present, and ice will build up along the cooling coils. Periodically defrost the ice (recommended when the ice level reaches a thickness of 1/4 to 1/2 inch). **NEVER USE A METAL INSTRUMENT** as you can damage the cooling coils, which will void your warranty and often cannot be repaired. To defrost, disconnect from the power supply, empty the contents, remove the drawers and provide a tray in the bottom to collect water. Use a hairdryer or other source of warm air to melt the ice (or leave the power off for 24 to 48 hours).

Appliance Care

Cleaning the Appliance:

NOTE: Use nonflammable cleaners.

Complete cleaning should be done at least once a year.

1. Unplug appliance or disconnect power.
2. Remove all stored items, then remove the drawers (see page 6).
3. Wash the inside walls and the drawers with a solution of mild detergent in warm water or 2 tbsp baking soda to 1 quart of warm water.
 - Do not use abrasive or harsh cleaners such as window sprays, scouring cleansers, flammable fluids, cleaning waxes, concentrated detergents, bleaches or cleansers containing petroleum products on plastic parts, interior and drawer liners or gaskets. Do not use paper towels, scouring pads, or other harsh cleaning tools. These can scratch or damage surfaces.
4. Rinse well and wipe dry with a clean soft cloth.
5. Use the appropriate cleaning method for the exterior finish.

Painted metal: Wash exteriors with a clean sponge or soft cloth and a mild detergent in warm water. Do not use abrasive or harsh cleaners, or cleaners designed for stainless steel. Dry thoroughly with a soft cloth. For additional protection against damage to painted metal exteriors, apply appliance wax (or auto paste wax) with a clean, soft cloth. Do not wax plastic parts.

Stainless steel (including interior of drawers): Routine cleaning can be accomplished by using warm water and a cloth. Dry with a towel or cloth to prevent water spots. Wipe in the directions of the brush lines. For cleaning that needs more power, use a mild detergent in warm water. Make sure to rinse the surface thoroughly to prevent staining and spotting. Towel dry to prevent water spots. As an alternative, use a commercial stainless steel cleaner, following directions carefully.

6. Replace drawers and plug in appliance or reconnect power.
7. Replace all stored items.
8. Every now and then wipe the condenser on the rear wall. Use a soft non-metal brush or vacuum cleaner.

Loss of Power:

After the power is restored, unplug the appliance for 30 minutes to allow the refrigeration system to equalize. During this time, keep the drawers closed as much as possible.

Remember: A full appliance stays cold longer than a partially filled one.

Moving the Appliance:

1. Remove stored items and tape any dividers in place.
2. Tape drawers shut.
3. Unplug the appliance.
4. Clean it thoroughly.
5. Tape the electrical cord to the cabinet.

When you get the appliance to its new location, refer to the *Installation* and *Temperature Selection* sections for information on installation and setting the controls.

Disposing of a Worn-Out Appliance:

- When an appliance wears out, dispose of it.
- Before disposing of your old refrigerator or freezer, take off the doors or remove the drawers so that children may not easily climb inside.
- The refrigerating system of the appliance is filled with refrigerant and insulating substances which should be separately treated and processed. Either have a licensed appliance repair company or dealer remove the appliance or call your local recycling office for the appropriate disposal.
- For the sake of environmental protection be careful not to damage the rear wall of the appliance (the condenser unit or the tubes when moving the freezer) or any part of the refrigerating system inside the appliance.

Troubleshooting

Try these solutions first to avoid the cost of an unnecessary service call.

The freezer will not operate:

- **Has a household fuse blown or circuit breaker tripped?**

Replace the fuse or reset the circuit breaker.

The compressor seems to run too much:

- **Is there excessive frost or a package preventing a drawer from closing completely?**

Defrost and clean the freezer, or move the package so the drawer closes properly.

- **Is the room temperature warmer than normal?**

The compressor will run longer under warm conditions. At normal room temperatures, it will run about 40% to 80% of the time. Under warmer conditions, it will run even more.

- **Have a large number of items just been added to the freezer?**

Adding a large amount warms the freezer. The compressor will normally run longer to cool the freezer back down.

- **Are the drawers closed completely?**

Push the drawers firmly shut. If they will not shut all the way, see *The drawers will not close completely* later in this section.

- **Are the condenser coils dirty (on some models)?**

This prevents heat transfer and makes the compressor work harder. Clean the condenser coils.

- **Are the drawer gaskets sealed all the way around?**

If not, contact a qualified technician.

Temperature is too warm:

- **Are the air vents blocked?**

This prevents cold air movement. Move objects away from the air vents.

- **Are the drawers opened often?**

Be aware that the freezer will warm when this occurs. In order to keep the freezer cool, try to get everything you need out of the freezer at once, keep items organized so they are easier to find, and close the drawer as soon as the items are removed.

The drawers will not close completely or are difficult to open:

- **Are packages blocking the drawer?**

Rearrange containers so that they fit more tightly and take up less space.

- **Are the drawers out of position?**

Reset the drawers back into their correct position (see page 6).

Limited Warranty

ONE YEAR LIMITED WARRANTY FOR APPLIANCE

Within the 48 contiguous United States, for one year from the date of purchase, when this appliance is operated and maintained according to instructions attached to or furnished with the product, warrantor will pay for factory-specified parts and repair labor to correct defects in materials or workmanship. Service must be provided by a designated service company. Outside the 48 states, all parts are warranted for one year from manufacturing defects. Plastic parts, shelves and cabinets are warranted to be manufactured to commercially acceptable standards, and are not covered from damage during handling or breakage.

5-YEAR COMPRESSOR WARRANTY

1. The compressor is covered for 5 years.
2. Replacement does not include labor.

ITEMS WARRANTOR WILL NOT PAY FOR:

1. Service calls to correct the installation of your appliance, to instruct you how to use your appliance, to replace or repair fuses or to correct wiring or plumbing.
2. Service calls to repair or replace appliance light bulbs or broken glass shelves. Consumable parts (such as filters) are excluded from warranty coverage.
3. Damage resulting from accident, alteration, misuse, abuse, fire, flood, acts of God, improper installation, installation not in accordance with electrical or plumbing codes, or use of products not approved by warrantor.
4. Replacement parts or repair labor costs for units operated outside the United States.
5. Repairs to parts or systems resulting from unauthorized modifications made to the appliance.
6. The removal and reinstallation of your appliance if it is installed in an inaccessible location or is not installed in accordance with published installation instructions.

DISCLAIMER OF IMPLIED WARRANTIES – LIMITATION OF REMEDIES

CUSTOMER'S SOLE AND EXCLUSIVE REMEDY UNDER THIS LIMITED WARRANTY SHALL BE PRODUCT REPAIR AS PROVIDED HEREIN. IMPLIED WARRANTIES, INCLUDING WARRANTIES OF MERCHANTABILITY OR FITNESS FOR A PARTICULAR PURPOSE, ARE LIMITED TO ONE YEAR. WARRANTOR SHALL NOT BE LIABLE FOR INCIDENTAL OR CONSEQUENTIAL DAMAGES. SOME STATES DO NOT ALLOW THE EXCLUSION OR LIMITATION OF INCIDENTAL OR CONSEQUENTIAL DAMAGES, OR LIMITATIONS ON THE DURATION OF IMPLIED WARRANTIES OF MERCHANTABILITY OR FITNESS, SO THESE EXCLUSIONS OR LIMITATIONS MAY NOT APPLY TO YOU. THIS WARRANTY GIVES YOU SPECIFIC LEGAL RIGHTS AND YOU MAY ALSO HAVE OTHER RIGHTS, WHICH VARY FROM STATE TO STATE.

WARNING: This product can expose you to chemicals including Nickel (Metallic) which is known to the State of California to cause cancer.

For more information go to www.P65Warnings.ca.gov

Note: Nickel is a component in all stainless steel and some other metal components.

Felix Storch, Inc.
770 Garrison Avenue
Bronx, New York 10474
Phone: (718) 893-3900
Fax: (844) 478-8799
www.summitappliance.com

For parts and accessory ordering,
troubleshooting and helpful hints,
visit: www.summitappliance.com/support

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